

Proteinase K Solution (Ready-To-Use)

REF / Cat. No.: ZY0007 15 ml
ZY0009 50 ml

Instructions for use

Intended use

Proteinase K Solution is a ready-to-use solution developed for enzymatic epitope retrieval on formalin-fixed tissue sections on slides. This procedure (sometimes called PIER, *Protease Induced Epitope Retrieval*) is primarily used in immunohistochemical staining procedures.

Proteinase K solution is intended for research use only.

Summary and explanations

Immunohistochemical staining procedures consist of sequential incubation steps with blocking solutions, antibodies and secondary reagents, enzymes and chromogenic substrates carried out on tissue sections. These tissue sections are mostly prepared out of formalin-fixed paraffin-embedded tissue blocks. Cellular structures are very effectively stabilised by formalin fixation which results in optimal morphological preservation of the sample.

On the other hand the formalin fixation leads to strong cross-links between proteins. This means that epitopes of antigens are being masked and often are no longer accessible for primary antibodies. In order to enable primary antibodies to bind to antigens the epitopes have to be recovered.

Enzymatic digestion with proteolytic enzymes (PIER) restores structures of the epitopes making them more accessible to specific antibodies. Heat induced epitope retrieval (HIER) in buffer solutions of different compositions and pH-values is another way of recovering epitopes. The primary antibody used determines the appropriate method.

Principle of the method

Proteinase K Solution is a ready-to-use Proteinase K solution in TBS buffer for enzymatic epitope retrieval.

Reagent provided

REF / Cat. No. ZY0007
15 ml Proteinase K Solution (Ready-To-Use)

REF / Cat. No. ZY0009
50 ml Proteinase K Solution (Ready-To-Use)

Storage and handling

The solution should be stored in aliquots at -20°C without further dilution. Under these conditions the solution is stable up to the expiry date indicated on the label. Do not use product after the expiry date.

A positive and a negative control have to be carried out in parallel to the test material. If you observe unusual staining or other deviations from the expected results which could possibly be caused by this reagent, please contact Zytomed Systems' technical support or your local distributor.

Precautions

Use by qualified personnel only. Wear protective clothing to avoid contact of the reagents with eye, skin or mucous membrane. In case of a reagent coming into contact with a sensitive area, wash the area with large amounts of water. Material safety data sheets (MSDS) are available upon request.

Reagent preparation

Proteinase K solution is ready-to-use and should be at room temperature prior to use.

Procedure

Proteinase K Solution is suitable for enzymatic epitope retrieval carried out after the dewaxing and rehydration of the tissue sections.

1. Apply 2-3 drops of ready-to-use Proteinase K Solution to deparaffinised and rehydrated tissue sections.
2. Incubate for 15 - 20 minutes at room temperature. The optimal incubation time needs to be elaborated by the operator.
3. Rinse carefully (3 x), first with distilled water followed by buffer.
4. Proceed with immunohistological staining as usual.

Quality control

We recommend carrying out a positive and a negative control with every staining run. The positive control permits the validation of appropriate processing of the sample. If the negative control has a positive result, this points to unspecific staining. Please refer to the instructions of the detection system for guidance on general quality control procedures.

Troubleshooting

If you observe unusual staining or other deviations from the expected results please read these instructions carefully, contact Zytomed Systems' technical support or your local distributor. Also refer to the instructions of the detection systems for guidance on general troubleshooting.

Expected results

During the reaction of the substrate with horse radish peroxidase or alkaline phosphatase in the presence of a chromogen, a coloured precipitate is formed at the location of the bound primary antibody. This reaction only takes place if the target antigen is existent in the tissue. The chromogen used determines the colour of the precipitate. The analysis is carried out using a light microscope.

Limitations of the procedure

Immunohistochemistry is a complex technique involving both histological and immunological detection methods. It requires a highly trained histotechnologist. Tissue processing and handling prior to immunostaining, for example variations in fixation and embedding or the inherent nature of the tissue can cause inconsistent results (Nadji and Morales, 1983). Inadequate counterstaining and mounting can influence the interpretation of the results. Zytomed Systems guarantees that the product will meet all requirements described from its shipping date until its expiry date, as long as the product is correctly stored and utilized. No additional guarantees can be given. Under no circumstances shall Zytomed System be liable for any damages arising out of the use of the reagent provided.

Performance characteristics

Zytomed Systems has conducted studies to evaluate the performance of the reagent. The product has been found to be suitable for the intended use.












Bibliography

Varma M, Linden MD, Amin MB Mod Pathol 12:472-478, 1999
 Nadji M and Morales AR Ann N.Y. Acad Sci 420:134-9, 1983

Thursday, 21 March 2013

ZY0007_9_Proteinase-K-A0313_english

Explanation of the symbols on the product label:

	Bestellnummer Catalog Number Reference du catalogue		Verwendbar bis Use By Utiliser jusque			Gebrauchsanweisung beachten Consult Instructions for use Consulter les instructions d'utilisation
	Chargenbezeichnung Batch Code Code du lot		Lagerungstemperatur Temperature Limitation Limites de température			Nur für Forschungszwecke For Research Use Only Pour la recherche uniquement
	In vitro Diagnostikum In Vitro Diagnostic Medical Device Dispositif médical de diagnostic in vitro		Achtung/Gefahr Warning/Danger Attention/Danger			Hersteller / Manufacturer / Fabricant Zytomed Systems GmbH Anhaltinerstraße 16 14163 Berlin, Germany Tel: (+49) 30-804 984 990 www.zytomed-systems.de
	Achtung/Gefahr Warning/Danger Attention/Danger		Achtung Warning Attention	