





Cap of the Pink PAP Pen:

Carefully pull off the cap in a straight line or with a slight twisting movement clockwise.

Tip of the Pink PAP Pen:

The tip of the Pink PAP Pen holds the fiber tip with which the water-repellent barrier is applied on the slide.

The tip is attached to the reservoir via a thread and must sit tightly closed on the reservoir. If necessary, please screw the tip tightly back onto the reservoir by turning it slightly clockwise.

Reservoir of the Pink PAP Pen:

The reservoir contains a water-repellent solution/reagent, which is applied to the slide via the fiber tip.

Shake vigorously or vortex before use, so that the reagent in the reservoir is mixed well, especially before the first use.

Fig. Design and handling of the Pink PAP Pen

**Limitations of the method**

Zytomed Systems guarantees that the product meets all requirements described from the time of dispatch to consumption, provided it is correctly stored and used. No further guarantees are given.

November 16, 2020

Rev: A1120

Doc: DBE\_ZUC064\_ZUC065

Explanations of the symbols on the product label:

<b>REF</b>	Bestellnummer Catalog Number Reference du catalogue		Verwendbar bis Use By Utiliser jusque		Gebrauchsanweisung beachten Consult Instructions for use Consulter les instructions d'utilisation
<b>LOT</b>	Chargenbezeichnung Batch Code Code du lot		Lagerungstemperatur Temperature Limitation Limites de température	<b>RUO</b>	Nur für Forschungszwecke For Research Use Only Pour la recherche uniquement
<b>IVD</b>	In vitro Diagnostikum In Vitro Diagnostic Medical Device Dispositif médical de diagnostic in vitro		Achtung/Gefahr Warning/Danger Attention/Danger		Hersteller / Manufacturer / Fabricant Zytomed Systems GmbH Anhaltinerstraße 16 14163 Berlin, Germany Tel: (+49) 30-804 984 990 <a href="http://www.zytomed-systems.de">www.zytomed-systems.de</a>
	Achtung/Gefahr Warning/Danger Attention/Danger		Achtung Warning Attention		